

Clean the Sash



For Non-motorized Sash:

- Remove armrest.
- Lower the sash until it is in the lowest position and creates a 10 cm (~4") opening between the sash and the diffuser.
- Open the front panel.
- Slide your hand through the opening and clean the inner side of the sash using cloth and IPA 70%.
- After cleaning, raise sash until it is in the sash ready position.
- Close the front panel.
- Reattach armrest.

For Motorized Sash:

- Remove armrest.
- Press MENU button, select SET MODE then choose MAINTENANCE MODE - this mode allows you to lower the sash further than usual.
- Lower the sash by pressing the DOWN button continuously until the sash is in the lowest position and creates a 10 cm (~4") opening between the sash and the diffuser.
- Open the front panel.
- Slide your hand through the opening and clean the inner side of the sash using cloth and IPA 70%.
- After cleaning, press the UP button continuously until the sash is in the sash ready position.
- Close the front panel.
- Reattach armrest.
- Set the cabinet back to NORMAL MODE

Product Lines	Size	Electrical
Labculture Class II, Type A2	LA2-3, LA2-4, LA2-5, LA2-6, LA2-4G1, LA2-6G1	A1/A2/A3/J7
Labculture Class II Low Noise	LA2-3L, LA2-4L, LA2-5L, LA2-6L	L1/L3
Labculture Class II, Type B2	LB2-3, LB2-4, LB2-5, LB2-6	B1/B2/B3/Y7
Cytoculture	CYT-4, CYT-6	A1/A2/A3
Airstream Class II (E-Series)	AC2-2E, AC2-3E, AC2-4E, AC2-5E, AC2-6E	E1/E2/E3
Airstream Class II (S-Series)	AC2-3S, AC2-4S, AC2-5S, AC2-6S	S1/S2/S3
Airstream DUO Class II (D-Series)	AC2-4D, AC2-6D	D1/D3
Airstream Class II (G-Series)	AC2-4G, AC2-6G	G1/G3
Airstream Class II, Type B2	AB2-3S, AB2-4S, AB2-5S, AB2-6S	S1/S2/S3
Airstream Max Class II	AC2-4L, AC2-6L	L1
Viva Universal Animal	VA2-4A, VA2-6A	A1/A2/A3
Labculture Plus Class II	LP2-4S, LP2-6S	S1
Labculture Plus Class II (D-Series)	LP2-4D	D1
Infinity Class II	FC2-4A, FC2-5A, FC2-6A	A1
Infinity Class II (Australia)	FC2-4K, FC2-6K	K1
Streamline Class II	SC2-4A, SC2-6A	A1/A2